



2024-12-17

Veritas Corporation
Sumitomo Hamamatsucho Bldg. 6th Floor
1-18-16 Hamamatsucho,
Minato-ku Tokyo 105-0013, JAPAN

To whom it may concern,

I am writing to inform you that, as part of continual improvements to our products, we are updating the recommended in-use storage duration for STEMdiff™ Hematopoietic Basal Medium (Catalog #05311). You are being notified because you have previously purchased a kit containing this product.

The new recommendation is as follows: If STEMdiff™ Hematopoietic Basal Medium is thawed but not used immediately, the medium may either be stored at 2-8°C for up to 4 weeks, or aliquoted and stored at -20°C ± 5°C. After thawing the aliquots, use immediately or store at 2-8°C for up to 2 weeks. Do not re-freeze the aliquots. Do not use beyond the shelf life of the basal medium.

The change in the in-use storage duration may apply to all batches of STEMdiff™ Hematopoietic Basal Medium (Catalog #05311). The Product Information Sheets have been updated to reflect this change. There is no impact on the product formulation, format, performance, or specifications as a result of this change.

Regards,



Brenda Clark
Manager, Quality Control, Stability